



K563XX ANHYDRIDE PURITY BATH

OPERATION AND INSTRUCTION MANUAL

REV A

Koehler Instrument Company, Inc.

1595 Sycamore Avenue • Bohemia, New York 11716-1796 • USA

Toll Free: 1-800-878-9070 (US only) • Tel: +1 631 589 3800 • Fax: +1 631 589 3815

<http://www.koehlerinstrument.com> • e-mail: info@koehlerinstrument.com

Petroleum Testing & Analysis Instrumentation • Custom Design & Manufacturing

CERTIFICATE OF CONFORMANCE

Anhydride Purity Bath

This certificate verifies that part number K563XX, Anhydride Purity Bath, was manufactured in conformance with the applicable standards set forth in this certification.

Specifications: **ASTM D3366**

This unit is tested before it leaves the factory, to ensure total functionality and compliance to the above specifications and ASTM standards. Test and inspection records are on file for verification.

June Kelly

Jesse Kelly
Application Engineer
Koehler Instrument Company

Table of Contents

1	INTRODUCTION.....	2
1.1	KOEHLER’S COMMITMENT TO OUR CUSTOMERS	2
1.2	RECOMMENDED RESOURCES AND PUBLICATIONS.....	2
1.3	INSTRUMENT SPECIFICATIONS	2
2	SAFETY INFORMATION AND WARNINGS	2
3	GETTING STARTED	3
3.1	PACKING LIST	3
3.2	UNPACKING	3
4	DESCRIPTIONS	3
4.1	TEMPERATURE CONTROLLER OPERATION	3
5	OPERATION.....	4
6	WIRING DIAGRAMS	5
7	REPLACEMENT PARTS.....	6
8	SERVICE	6
9	STORAGE	6
10	WARRANTY	6
11	RETURNED GOODS POLICY.....	6
	NOTES	7
	NOTES	8

1 Introduction

The K563XX Anhydride Purity Bath is an electrically heated aluminum block bath with a microprocessor-based temperature controller. Molten samples of maleic or phthalic anhydride are compared to Platinum-Cobalt color standards to determine the sample purity and the qualitative stability in the presence of contaminants.

The aluminum block bath is equipped with over temperature protection circuitry and a dual LED temperature display. The heating unit provides temperature stability, heating rates, and minimal temperature gradients which exceed ASTM specifications, and is housed in an insulated steel cabinet with a chemically-resistant painted finish. Up to six samples can be tested at one time using Nessler Tubes.

This manual provides important information regarding safety, technical reference, installation requirements, operating condition specifications, user facility resource requirements, and operating instructions for the Cloud and Pour Point Bath. This manual should also be used in conjunction with applicable published laboratory procedures. Information on these procedures is given in section 1.2.

1.1 Koehler's Commitment to Our Customers

Providing quality testing instrumentation and technical support services for research and testing laboratories has been our specialty for more than 50 years. At Koehler, the primary focus of our business is providing you with the full support of your laboratory testing needs. Our products are backed by our staff of technically knowledgeable, trained specialists who are experienced in both petroleum products testing and instrument service to better understand your requirements and provide you with the best solutions. You can depend on Koehler for a full range of accurate and reliable instrumentation as well as support for your laboratory testing programs. Please do not hesitate to contact us at any time with your inquiries about equipment, tests, or technical support.

Toll Free: 1-800-878-9070 (US only)
Tel: +1 631 589 3800
Fax: +1 631 589 3815

Email: info@koehlerinstrument.com
<http://www.koehlerinstrument.com>

1.2 Recommended Resources and Publications

1. American Society for Testing and Materials (ASTM)
100 Barr Harbor Drive
West Conshohocken, Pennsylvania 19428-2959, USA
Tel: +1 610 832 9500
Fax: +1 610 832 9555
<http://www.astm.org>
email: service@astm.org

ASTM Publication:

- ASTM D3366: Color of Maleic Anhydride and Phthalic Anhydride in the Molten State and after Heating (Platinum-Cobalt Scale)

1.3 Instrument Specifications

Model Number	Voltage	Frequency
K56300	115V	60Hz
K56390	220-240V	50Hz/60Hz

2 Safety Information and Warnings

Safety Considerations. The use of this equipment may involve *hazardous* materials and operations. This manual does not purport to address all of the safety problems associated with the use of this equipment. It is the responsibility of any user of this equipment to investigate, research, and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

Equipment Modifications and Replacement Parts. Any modification or alteration of this equipment from that of factory specifications is not recommended voids the manufacturer warranty, product safety, performance specifications, and/or certifications whether specified or implied, and may result in personal injury and/or property loss. Replacement parts must be O.E.M. exact replacement equipment.

Unit Design. This equipment is specifically designed for use in accordance with the applicable standard test methods listed in section 1.2 of this manual. The use of this

equipment in accordance with any other test procedures, or for any other purpose, is not recommended and may be extremely hazardous.

This unit is NOT explosion-proof and proper precautionary measures should be taken during operation. Avoid using this unit in an environment with flammable or explosive vapors.

Chemical Reagents Information. Chemicals and reagents used in performing the test may exhibit potential hazards. Any user must be familiarized with the possible dangers before use. We also recommend consulting the Material Data and Safety Sheet (MSDS) on each chemical reagent for additional information. MSDS information can be easily located on the internet at <http://siri.uvm.edu> or <http://www.sigma-aldrich.com>.

3 Getting Started

The instructions for preparing the equipment assume that the user is aware of the contents of this document, which lists the warranty conditions and important precautions.

3.1 Packing List

- K563XX Anhydride Purity Bath
- K563XX-Manual Anhydride Purity Bath Operation and Instruction Manual

Accessories (To be purchased separately)

- K56306 Nessler Tubes (set of 6)

3.2 Unpacking

This unit comes packed in a crate. Remove the bath from the crate and place on a firm, level table/

Unwrap the unit and wipe the shell with a soft cloth to remove smudges caused by wrapping material.

Connect the line cord to a properly fused and grounded receptacle with the correct voltage as indicated in section 1.3.

WARNING: For safety, disconnect the power when performing any maintenance and/or cleaning. Do **NOT** turn the power on unless the

bath is filled with the proper medium; otherwise, damage may occur to the unit and the warranty will be void.

4 Descriptions

4.1 Temperature Controller Operation

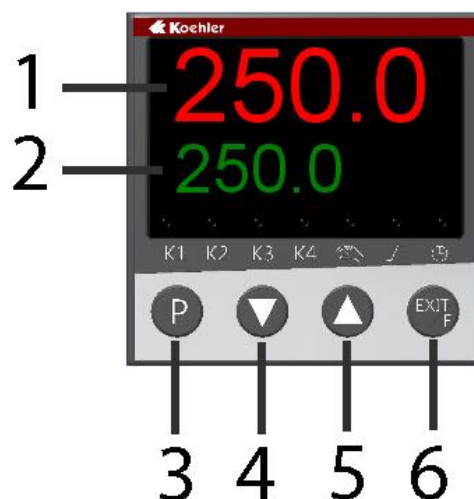


Figure 4. Temperature Controller

- 1. Process Temperature Display.** The upper red LED display shows the process temperature as read from the RTD probe.
- 2. Set Point Temperature Display.** The lower green LED display shows the set point temperature of the controller.
- 3. Programming Key.** Permits scrolling through controller menu parameters. One Level Forward
- 4. Down Key.** Used to decrease the set point temperature and to decrease or change parameter values when programming the temperature controller.
- 5. Up Key.** Used to increase the set point temperature and to increase or change parameter values when programming the temperature controller.
- 6. Exit / Function Key.** This key is used to exit or leave a level. One level backward

IMPORTANT NOTE: The digital temperature controller for the unit comes pre-programmed from the Koehler factory. Please do NOT attempt to re-program the digital temperature controller as this will void the product warranty. If assistance is required, please do not hesitate to contact the Koehler technical service department.

Setting the Temperature. Set the desired operating temperature by adjusting the set point with the up and down keys. The set point will be displayed in the lower green Set Point LED display and the actual temperature will be displayed in the upper red Process LED display. Please allow the instrument to fully equilibrate before proceeding with any testing.

Temperature Calibration. This routine allows the digital temperature controller to be calibrated to a certified thermometer.

- a. Use a certified calibrated measuring device to acquire the temperature. Calculate the difference between the measuring device and the Process value displayed on the controller.
- b. Press the program key two times until **PCT** is displayed in the lower green LED display. Press the DOWN key. CAL will display on the lower green display. If there is a value observed in the upper red LED display, add it to the calculated difference obtained in the previous step. This is the offset value.
- c. Press the Program Key. The lower green display will flash. Use the up or down keys to adjust to the new calibration offset value on the upper red display calculated in the previous step. When the value has been entered, the controller will automatically store the value. The lower green display will stop flashing. If further adjustments are necessary, press the Program Key again. Resume regular operations by pressing the Exit / Function key two times. Verify if the new calibration is correct by observing the upper red display and comparing the value with the calibrated reference device.

Auto Tune. This routine allows the digital temperature control to learn the heating parameters needed for any particular set point temperature. This operation should be done when installing a new unit, after replacing or changing the bath medium type, or utilizing a different temperature set point 20% different from the previously used set point temperature.

- a. Set the operating temperature to the desired setting.
- b. Press the up and down arrow buttons simultaneously for about 5 seconds. When Auto Tune is active, the lower green LED display will blink **TUNE**. Auto Tune will automatically toggle off when the set point temperature is reached. Auto tune can be terminated by pressing the up & down buttons simultaneously again.


5 Operation

Be sure to read the safety and hazard warnings, the installation procedure and any of the standard test methods mentioned in the Introduction of this manual before operating this unit.

Turn ON the line switch. This switch is located on the bottom right-hand corner of the unit.

Insert a calibrated thermometer into the aluminum block thermometer port. Observe the display panel of the digital temperature control.

6 Wiring Diagrams

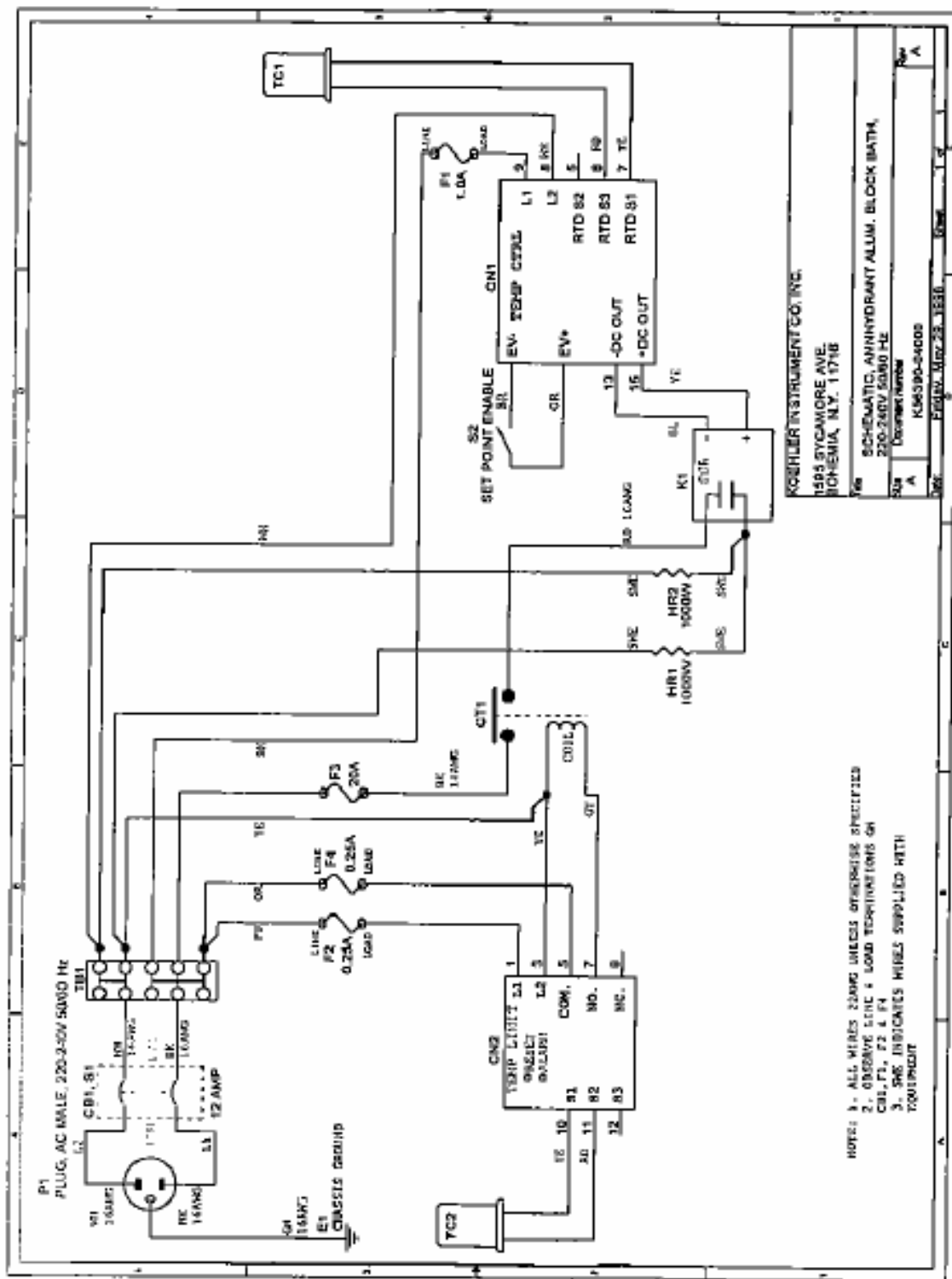

7 Replacement Parts

Part Number	Description
275-103-044	Electronic Temperature Controller
220-240-014	Cartridge Heater, 1000W 240V
265—203-001	Temperature Probe, Type K

8 Service

Under normal operating conditions and with routine maintenance, the K563XX Anhydride Purity Bath should not require service. Any service problem can be quickly resolved by contacting Koehler's technical service department either by letter, phone, fax, or email. In order to assure the fastest possible service, please provide us with the following information.

Model Number: _____

Serial Number: _____

Date of Shipment: _____

9 Storage

This laboratory test instrument is equipped with electrical components. Storage facilities should be consistent with an indoor laboratory environment. This testing equipment should not be subjected to extremes of temperature and/or moisture.

This equipment was shipped from the factory in a corrugated cardboard container. If long term storage is anticipated, re-packing the instrument in a water-resistant container is recommended to ensure equipment safety and longevity.


10 Warranty

We, at Koehler, would like to thank you for your equipment purchase, which is protected by the following warranty. If within one (1) year from the date of receipt, but no longer than fifteen (15) months from the date of shipment, Koehler equipment fails to perform properly because of defects in materials or workmanship, Koehler Instrument Company, Inc. will repair or, at its sole discretion, replace the equipment without charge F.O.B. its plant, provided the equipment has been properly installed, operated, and maintained. Koehler Instrument Company must

be advised in writing of the malfunction and authorize the return of the product to the factory. The sole responsibility of Koehler Instrument Company and the purchaser's exclusive remedy for any claim arising out of the purchase of any product is the repair or replacement of the product. In no event shall the cost of the purchaser's remedy exceed the purchase price, nor shall Koehler Instrument Company be liable for any special, indirect, incidental, consequential, or exemplary damages. KOEHLER INSTRUMENT COMPANY, INC. DISCLAIMS ALL OTHER WARRANTIES, EXPRESSED OR IMPLIED, INCLUDING ANY IMPLIED WARRANTIES OF FITNESS FOR A PARTICULAR PURPOSE. Please save the shipping carton in the event the equipment needs to be returned to the factory for warranty repair. If the carton is discarded, it will be the purchaser's responsibility to provide an appropriate shipping carton.

11 Returned Goods Policy

To return products for credit or replacement, please contact Koehler Customer Service with your purchase order number, our packing list/invoice number, the item(s) to be returned and the reason for the return. You will be issued a Returned Authorization (RA) number, which must be prominently displayed on the shipping container when you return the material to our plant. Shipping containers without an RA number prominently displayed with will be returned to the sender. Goods must be returned freight prepaid. Returns will be subject to a restocking charge, the application of which will depend upon the circumstances necessitating the return. Some returns cannot be authorized, including certain products purchased from outside vendors for the convenience of the customer, products manufactured on special order, products shipped from the factory past ninety (90) days, and products which have been used or modified in such a way that they cannot be returned to stock for future sale.

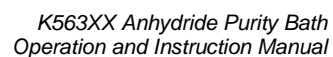[illegible]



*K563XX Anhydride Purity Bath
Operation and Instruction Manual*